

Device Tools Windows Help

- Connect..
- Disconnect
- Exit

Connect to the instrument by going to 'Device' and 'Connect'

Find your instrument's IP address and click 'Connect'

Connect

Network Connection Serial Communication Direct Hardware Access

**Active Remote Connection:**

Name	IP Address
800N2682	192.168.0.36

Connect Cancel

Name	Size	Type	Date Modified
------	------	------	---------------

- 800N2682
  - Application Data
  - ARCHIVE
  - BINARIES
  - BRUKER
  - LBOOT
  - My Documents
  - Network
  - Program Files
  - Recycled
  - Temp

A file explorer window will open up once you are connected

- My Computer
  - PRL (A:)
  - (C:)
  - \\Mac\Google Drive\ (V:)
  - \\Mac\Dropbox\ (W:)
  - \\Mac\Photo Library\ (X:)
  - \\Mac\iCloud\ (Y:)
  - \\Mac\Home\ (Z:)

\BRUKER

- 800N2682
  - Application Data
  - ARCHIVE
  - BINARIES
  - BRUKER**
    - Autold
    - Data
    - Documents and Settings
    - EDX Databases
    - lost+found
    - package
    - Setup
    - System
    - TitanCals
  - LBOOT
  - My Documents
  - Network
  - Program Files
  - Recycled
  - Temp

Name	Size	Type	Date Modified
Autold		File Fold...	1/1/06 2:00 PM
Data		File Fold...	10/21/14 11:35 AM
Documents and Settings		File Fold...	1/1/06 6:00 AM
EDX Databases		File Fold...	1/1/06 2:00 PM
lost+found		File Fold...	1/11/06 12:59 PM
package		File Fold...	1/1/06 2:00 PM
BruckerInstallerLog0.txt	141 KB	Text Do...	12/6/16 3:04 AM
Cal DualMining Context ...	8 KB	SNPSH...	5/30/16 8:10 AM
Cargill.aen	5 KB	AEN File	6/20/16 11:08 AM
eventlog0.txt	1,505 KB	Text Do...	1/1/06 2:01 PM
eventlog1.txt	721 KB	Text Do...	1/11/06 9:11 PM
eventlog2.txt	931 KB	Text Do...	1/1/06 9:01 PM
GeoChem.aen	22 KB	AEN File	4/11/16 10:21 AM
InstrumentDef.idf	34 KB	IDF File	1/11/06 9:00 PM
MudrockMajor.aen	5 KB	AEN File	7/13/16 1:53 PM
Mudrock-Majors.aen	5 KB	AEN File	5/5/16 1:16 PM
MudrockTrace.aen	5 KB	AEN File	7/13/16 3:18 PM
Mudrock-Trace.aen	5 KB	AEN File	5/5/16 3:01 PM
Obsidian.aen	5 KB	AEN File	7/13/16 3:27 PM
plantCal.aen	5 KB	AEN File	12/6/16 2:57 AM
PreciousMetals.aen	19 KB	AEN File	11/2/15 2:04 PM
RoHS.alt	47 KB	ALT File	11/2/15 2:04 PM
SmartGrade.aen	52 KB	AEN File	4/11/16 10:21 AM
Soil.aen	21 KB	AEN File	4/11/16 10:21 AM
SOILTEST50.aen	5 KB	AEN File	4/30/16 4:18 PM
StdAlloy.aen	1,411 KB	AEN File	4/11/16 10:22 AM
tracelog0.txt	523 KB	Text Do...	1/1/06 9:01 PM

Go to the BRUKER folder

- My Computer
  - PRL (A:)
  - (C:)
  - \\Mac\Google Drive\ (V:)
  - \\Mac\Dropbox\ (W:)
  - \\Mac\Photo Library\ (X:)
  - \\Mac\iCloud\ (Y:)
  - \\Mac\Home\ (Z:)



\BRUKER

Name	Size	Type	Date Modified
Autold		File Fold...	1/1/06 2:00 PM
Data		File Fold...	10/21/14 11:35 AM
Documents and Settings		File Fold...	1/1/06 6:00 AM
EDX Databases		File Fold...	1/1/06 2:00 PM
lost+found		File Fold...	1/11/06 12:59 PM
package		File Fold...	1/1/06 2:00 PM
Setup		File Fold...	1/1/06 2:00 PM
System		File Fold...	1/1/06 2:00 PM
TitanCals		File Fold...	10/27/14 12:26 PM
BrukerInstaller.xls	1 KB	Microsof...	10/21/14 11:35 AM
BrukerInstallerLog0.txt	141 KB	Text Do...	12/6/16 3:04 AM
Cal DualMining Context ...	8 KB	SNPSH...	5/30/16 8:10 AM
Cargill.aen	5 KB	AEN File	6/20/16 11:08 AM
eventlog0.txt	1,505 KB	Text Do...	1/1/06 2:01 PM
eventlog1.txt	721 KB	Text Do...	1/11/06 9:11 PM
eventlog2.txt	931 KB	Text Do...	1/1/06 9:01 PM
GeoChem.aen	22 KB	AEN File	4/11/16 10:21 AM
InstrumentDef.idf	34 KB	IDF File	1/11/06 9:00 PM
MudrockMajor.aen	5 KB	AEN File	7/13/16 1:53 PM
Mudrock-Majors.aen	5 KB	AEN File	5/5/16 1:16 PM
MudrockTrace.aen	5 KB	AEN File	7/13/16 3:18 PM
Mudrock-Trace.aen	5 KB	AEN File	5/5/16 3:01 PM
Obsidian.aen	5 KB	AEN File	7/13/16 3:27 PM
plantCal.aen	5 KB	AEN File	12/6/16 2:57 AM
PreciousMetals.aen	19 KB	AEN File	11/2/15 2:04 PM
RoHS.alt	47 KB	ALT File	11/2/15 2:04 PM
SmartGrade.aen	52 KB	AEN File	4/11/16 10:21 AM
Soil.aen	21 KB	AEN File	4/11/16 10:21 AM
SOILTEST50.aen	5 KB	AEN File	4/30/16 4:18 PM
StdAlloy.aen	1,411 KB	AEN File	4/11/16 10:22 AM
tracelog0.txt	523 KB	Text Do...	1/1/06 9:01 PM

The contents of your instrument will be here, including calibrations and data

My Computer

- ▶ PRL (A:)
- ▶ (C:)
- ▶ \\Mac\Google Drive\ (V:)
- ▶ \\Mac\Dropbox\ (W:)
- ▶ \\Mac\Photo Library\ (X:)
- ▶ \\Mac\iCloud\ (Y:)
- ▶ \\Mac\Home\ (Z:)

\BRUKER\Data

- 800N2682
  - Application Data
  - ARCHIVE
  - BINARIES
  - BRUKER
    - Autold
    - Data
    - Documents and Settings
    - EDX Databases
    - lost+found
    - package
    - Setup
    - System
    - TitanCals
  - LBOOT
  - My Documents
  - Network
  - Program Files
  - Recycled
  - Temp

Name	Size	Type	Date Modified
00010-Mudrock-Majors...	8 KB	PDZ File	6/6/16 2:54 AM
00012-Cargill.pdz	8 KB	PDZ File	6/6/16 2:54 AM
00025-Mudrock-Majors...	16 KB	PDZ File	6/13/16 6:43 PM
00026-Mudrock-Trace.p...	16 KB	PDZ File	6/13/16 6:43 PM
00039-Mudrock-Trace.p...	16 KB	PDZ File	5/28/16 5:15 AM
00040-Mudrock-Trace.p...	16 KB	PDZ File	5/28/16 5:16 AM
00048-Mudrock-Majors...	16 KB	PDZ File	7/8/16 1:51 PM
00052-Spectrometer Mo...	17 KB	PDZ File	7/9/16 8:23 PM
00054-Spectrum Only.pdz	8 KB	PDZ File	5/3/16 7:25 PM
00061-Mudrock-Majors...	8 KB	PDZ File	7/18/16 7:45 AM
00063-MudrockMajor.pdz	8 KB	PDZ File	7/18/16 7:47 AM
00064-MudrockMajor.pdz	8 KB	PDZ File	7/18/16 7:49 AM
00066-MudrockMajor.pdz	8 KB	PDZ File	7/18/16 7:52 AM
00070-MudrockMajor.pdz	8 KB	PDZ File	7/18/16 7:55 AM
00071-MudrockMajor.pdz	8 KB	PDZ File	7/18/16 7:55 AM
00071-Mudrock-Trace.p...	8 KB	PDZ File	6/5/16 1:26 AM
00072-MudrockMajor.pdz	8 KB	PDZ File	7/18/16 7:56 AM
00075-MudrockMajor.pdz	8 KB	PDZ File	7/18/16 7:58 AM
00076-MudrockMajor.pdz	8 KB	PDZ File	7/18/16 7:58 AM
00079-MudrockMajor.pdz	8 KB	PDZ File	7/18/16 8:02 AM
00080-MudrockMajor.pdz	8 KB	PDZ File	7/18/16 8:03 AM
00081-MudrockMajor.pdz	8 KB	PDZ File	7/18/16 8:04 AM
00082-Spectrum Only.pdz	8 KB	PDZ File	5/4/16 2:07 PM
00083-Spectrum Only.pdz	8 KB	PDZ File	5/4/16 2:08 PM
00084-MudrockMajor.pdz	8 KB	PDZ File	7/18/16 8:07 AM
00085-MudrockMajor.pdz	8 KB	PDZ File	7/18/16 8:10 AM
00087-Spectrum Only.pdz	8 KB	PDZ File	5/4/16 2:14 PM
00088-MudrockMajor.pdz	8 KB	PDZ File	7/18/16 8:11 AM
00089-MudrockMajor.pdz	8 KB	PDZ File	7/18/16 8:11 AM

Navigate to the Data folder

- My Computer
  - PRL (A:)
  - (C:)
  - \\Mac\Google Drive\ (V:)
  - \\Mac\Dropbox\ (W:)
  - \\Mac\Photo Library\ (X:)
  - \\Mac\iCloud\ (Y:)
  - \\Mac\Home\ (Z:)



\BRUKER\Data

Name	Size	Type	Date Modified
800N2682			
Application Data			
ARCHIVE			
BINARIES			
BRUKER			
Autold			
Data			
Documents and Settings			
EDX Databases			
lost+found			
package			
Setup			
System			
TitanCals			
LBOOT			
My Documents			
Network			
Program Files			
Recycled			
Temp			
00010-Mudrock-Majors...	8 KB	PDZ File	6/6/16 2:54 AM
00012-Cargill.pdz	8 KB	PDZ File	6/6/16 2:54 AM
00025-Mudrock-Majors...	16 KB	PDZ File	6/13/16 6:43 PM
00026-Mudrock-Trace.p...	16 KB	PDZ File	6/13/16 6:43 PM
00039-Mudrock-Trace.p...	16 KB	PDZ File	5/28/16 5:15 AM
00040-Mudrock-Trace.p...	16 KB	PDZ File	5/28/16 5:16 AM
00040-Spectrum Only.pdz	16 KB	PDZ File	7/8/16 4:54 PM
00047-GeoChem.pdz	9 KB	PDZ File	5/3/16 6:54 PM
00047-Spectrometer Mo...	17 KB	PDZ File	7/9/16 8:21 PM
00051-Spectrometer Mo...	17 KB	PDZ File	7/9/16 8:23 PM
00052-Spectrometer Mo...	17 KB	PDZ File	7/9/16 8:23 PM
00054-Spectrum Only.pdz	8 KB	PDZ File	5/3/16 7:25 PM
00061-Mudrock-Majors...	8 KB	PDZ File	7/18/16 7:45 AM
00063-MudrockMajor.pdz	8 KB	PDZ File	7/18/16 7:47 AM
00064-MudrockMajor.pdz	8 KB	PDZ File	7/18/16 7:49 AM
00066-MudrockMajor.pdz	8 KB	PDZ File	7/18/16 7:52 AM
00070-MudrockMajor.pdz	8 KB	PDZ File	7/18/16 7:55 AM
00071-MudrockMajor.pdz	8 KB	PDZ File	7/18/16 7:55 AM
00071-Mudrock-Trace.p...	8 KB	PDZ File	6/5/16 1:26 AM
00072-MudrockMajor.pdz	8 KB	PDZ File	7/18/16 7:56 AM
00075-MudrockMajor.pdz	8 KB	PDZ File	7/18/16 7:58 AM
00076-MudrockMajor.pdz	8 KB	PDZ File	7/18/16 7:58 AM
00079-MudrockMajor.pdz	8 KB	PDZ File	7/18/16 8:02 AM
00080-MudrockMajor.pdz	8 KB	PDZ File	7/18/16 8:03 AM
00081-MudrockMajor.pdz	8 KB	PDZ File	7/18/16 8:04 AM
00082-Spectrum Only.pdz	8 KB	PDZ File	5/4/16 2:07 PM
00083-Spectrum Only.pdz	8 KB	PDZ File	5/4/16 2:08 PM
00084-MudrockMajor.pdz	8 KB	PDZ File	7/18/16 8:07 AM
00085-MudrockMajor.pdz	8 KB	PDZ File	7/18/16 8:10 AM
00087-Spectrum Only.pdz	8 KB	PDZ File	5/4/16 2:14 PM
00088-MudrockMajor.pdz	8 KB	PDZ File	7/18/16 8:11 AM

In this folder will be all the spectra you have taken. If a USB is in your instrument, the data will be saved there instead

My Computer

- PRL (A:)
- (C:)
- \\Mac\Google Drive\ (V:)
- \\Mac\Dropbox\ (W:)
- \\Mac\Photo Library\ (X:)
- \\Mac\iCloud\ (Y:)
- \\Mac\Home\ (Z:)



\BRUKER\Data

Name	Size	Type	Date Modified
01130-MudrockTrace.pdz	9 KB	PDZ File	4/14/17 12:16 PM
01131-MudrockMajor.pdz	8 KB	PDZ File	4/14/17 12:17 PM
01132-MudrockMajor.pdz	8 KB	PDZ File	4/14/17 12:18 PM
01133-MudrockTrace.pdz	9 KB	PDZ File	4/14/17 12:21 PM
01134-MudrockMajor.pdz	8 KB	PDZ File	4/14/17 12:22 PM
01135-MudrockMajor.pdz	8 KB	PDZ File	4/14/17 12:23 PM
01136-MudrockTrace.pdz	9 KB	PDZ File	4/14/17 12:24 PM
01137-MudrockMajor.pdz	8 KB	PDZ File	4/14/17 12:26 PM
01138-Spectrum Only.pdz	8 KB	PDZ File	4/14/17 12:48 PM
01139-Spectrum Only.pdz	8 KB	PDZ File	4/14/17 12:49 PM
01140-MudrockMajor.pdz	8 KB	PDZ File	4/14/17 1:11 PM
01141-MudrockTrace.pdz	9 KB	PDZ File	4/14/17 1:12 PM
01142-MudrockMajor.pdz	8 KB	PDZ File	4/14/17 1:18 PM
01143-MudrockMajor.pdz	8 KB	PDZ File	4/14/17 1:34 PM
01144-MudrockMajor.pdz	8 KB	PDZ File	4/14/17 1:35 PM
01145-MudrockMajor.pdz	8 KB	PDZ File	4/14/17 1:36 PM
01146-MudrockMajor.pdz	8 KB	PDZ File	4/14/17 1:38 PM
01147-MudrockTrace.pdz	9 KB	PDZ File	4/14/17 1:39 PM
01148-MudrockTrace.pdz	9 KB	PDZ File	4/14/17 1:40 PM
01149-MudrockTrace.pdz	9 KB	PDZ File	4/14/17 1:42 PM
01150-MudrockTrace.pdz	9 KB	PDZ File	4/14/17 1:43 PM
01151-Spectrometer Mo...	8 KB	PDZ File	1/1/06 2:01 PM
01152-Spectrometer Mo...	8 KB	PDZ File	1/1/06 2:02 PM
01153-Spectrometer Mo...	8 KB	PDZ File	1/1/06 2:03 PM
01154-MudrockTrace.pdz	9 KB	PDZ File	1/1/06 9:01 PM
01155-MudrockTrace.pdz	9 KB	PDZ File	1/1/06 9:01 PM
01156-MudrockTrace.pdz	9 KB	PDZ File	1/1/06 9:02 PM
01157-MudrockTrace.pdz	9 KB	PDZ File	1/1/06 9:05 PM
01158-MudrockTrace.pdz	9 KB	PDZ File	1/1/06 9:06 PM
01159-MudrockTrace.pdz	9 KB	PDZ File	1/1/06 9:06 PM
Results.csv	520 KB	Microsof...	1/1/06 9:06 PM

This folder also has a 'Results.csv' file, which is a spreadsheet of your quantitative results

My Computer

- ▶ PRL (A:\)
- ▶ (C:\)
- ▶ \\Mac\Google Drive\ (V:\)
- ▶ \\Mac\Dropbox\ (W:\)
- ▶ \\Mac\Photo Library\ (X:\)
- ▶ \\Mac\iCloud\ (Y:\)
- ▶ \\Mac\Home\ (Z:\)

\BRUKER

- 800N2682
  - Application Data
  - ARCHIVE
  - BINARIES
  - BRUKER
    - Autold
    - Data
    - Documents and Settings
    - EDX Databases
    - lost+found
    - package
    - Setup
    - System
    - TitanCals
  - LBOOT
  - My Documents
  - Network
  - Program Files
  - Recycled
  - Temp

Name	Size	Type	Date Modified
Autold		File Fold...	1/1/06 2:00 PM
Data		File Fold...	10/21/14 11:35 AM
Documents and Settings		File Fold...	1/1/06 6:00 AM
EDX Databases		File Fold...	1/1/06 2:00 PM
lost+found		File Fold...	1/11/06 12:59 PM
package		File Fold...	1/1/06 2:00 PM
Setup		File Fold...	1/1/06 2:00 PM
System		File Fold...	1/1/06 2:00 PM
TitanCals		File Fold...	10/27/14 12:26 PM
BrukerInstaller.xls	1 KB	Microsof...	10/21/14 11:35 AM
BrukerInstallerLog0.txt	141 KB	Text Do...	12/6/16 3:04 AM
Cal DualMining Context ...	8 KB	SNPSH...	5/30/16 8:10 AM
Cargill.aen	5 KB	AEN File	6/20/16 11:08 AM
eventlog0.txt	1,505 KB	Text Do...	1/1/06 2:01 PM
eventlog1.txt	721 KB	Text Do...	1/11/06 9:11 PM
eventlog2.txt	931 KB	Text Do...	1/1/06 9:01 PM
GeoChem.aen	22 KB	AEN File	4/11/16 10:21 AM
InstrumentDef.idf	34 KB	IDF File	1/11/06 9:00 PM
MudrockMajor.aen	5 KB	AEN File	7/13/16 1:53 PM
Mudrock-Majors.aen	5 KB	AEN File	5/5/16 1:16 PM
MudrockTrace.aen	5 KB	AEN File	7/13/16 3:18 PM
Mudrock-Trace.aen	5 KB	AEN File	5/5/16 3:01 PM
Obsidian.aen	5 KB	AEN File	7/13/16 3:27 PM
plantCal.aen	5 KB	AEN File	12/6/16 2:57 AM
PreciousMetals.aen	19 KB	AEN File	11/2/15 2:04 PM
RoHS.alt	47 KB	ALT File	11/2/15 2:04 PM
SmartGrade.aen	52 KB	AEN File	4/11/16 10:21 AM
Soil.aen	21 KB	AEN File	4/11/16 10:21 AM
SOILTEST50.aen	5 KB	AEN File	4/30/16 4:18 PM
StdAlloy.aen	1,411 KB	AEN File	4/11/16 10:22 AM
tracelog0.txt	523 KB	Text Do...	1/1/06 9:01 PM

Go back one level, and select the 'Data' folder

- My Computer
  - PRL (A:)
  - (C:)
  - \\Mac\Google Drive\ (V:)
  - \\Mac\Dropbox\ (W:)
  - \\Mac\Photo Library\ (X:)
  - \\Mac\iCloud\ (Y:)
  - \\Mac\Home\ (Z:)



\BRUKER

- 800N2682
  - Application Data
  - ARCHIVE
  - BINARIES
  - BRUKER
    - Autold
    - Data
    - Documents and Settings
    - EDX Databases
    - lost+found
    - package
    - Setup
    - System
    - TitanCals
  - LBOOT
  - My Documents
  - Network
  - Program Files
  - Recycled
  - Temp

Name	Size	Type	Date Modified
Autold		File Fold...	1/1/06 2:00 PM
Data		File Fold...	10/21/14 11:35 AM
Documents and Settings		File Fold...	1/1/06 6:00 AM
EDX Databases		File Fold...	1/1/06 2:00 PM
lost+found			
package			
Setup			
System			
TitanCals			
BrukerInstallerLogo.txt	141 KB	Text Do...	12/6/16 3:04 AM
Cal DualMining Context ...	8 KB	SNPSH...	5/30/16 8:10 AM
Cargill.aen	5 KB	AEN File	6/20/16 11:08 AM
eventlog0.txt	1,505 KB	Text Do...	1/1/06 2:01 PM
eventlog1.txt	721 KB	Text Do...	1/11/06 9:11 PM
eventlog2.txt	931 KB	Text Do...	1/1/06 9:01 PM
GeoChem.aen	22 KB	AEN File	4/11/16 10:21 AM
InstrumentDef.idf	34 KB	IDF File	1/11/06 9:00 PM
MudrockMajor.aen	5 KB	AEN File	7/13/16 1:53 PM
Mudrock-Majors.aen	5 KB	AEN File	5/5/16 1:16 PM
MudrockTrace.aen	5 KB	AEN File	7/13/16 3:18 PM
Mudrock-Trace.aen	5 KB	AEN File	5/5/16 3:01 PM
Obsidian.aen	5 KB	AEN File	7/13/16 3:27 PM
plantCal.aen	5 KB	AEN File	12/6/16 2:57 AM
PreciousMetals.aen	19 KB	AEN File	11/2/15 2:04 PM
RoHS.alt	47 KB	ALT File	11/2/15 2:04 PM
SmartGrade.aen	52 KB	AEN File	4/11/16 10:21 AM
Soil.aen	21 KB	AEN File	4/11/16 10:21 AM
SOILTEST50.aen	5 KB	AEN File	4/30/16 4:18 PM
StdAlloy.aen	1,411 KB	AEN File	4/11/16 10:22 AM
tracelog0.txt	523 KB	Text Do...	1/1/06 9:01 PM

With 'Data' selected, click 'Download'

- My Computer
  - PRL (A:)
  - (C:)
  - \\Mac\Google Drive\ (V:)
  - \\Mac\Dropbox\ (W:)
  - \\Mac\Photo Library\ (X:)
  - \\Mac\iCloud\ (Y:)
  - \\Mac\Home\ (Z:)

BRUKER

- 800N2682
  - Application Data
  - ARCHIVE
  - BINARIES
  - BRUKER
    - Autold
    - Data
    - Documents and Settings
    - EDX Databases
    - lost+found
    - package
    - Setup
    - System
    - TitanCals
  - LBOOT
  - My Documents
  - Network
  - Program Files
  - Recycled
  - Temp

Name	Size	Type	Date Modified
Autold		File Fold...	1/1/06 2:00 PM
Data		File Fold...	10/21/14 11:35 AM
Documents and Settings		File Fold...	1/1/06 6:00 AM
EDX Databases		File Fold...	1/1/06 2:00 PM
lost+found		File Fold...	1/11/06 12:59 PM
package		File Fold...	1/1/06 2:00 PM
Setup		File Fold...	1/1/06 2:00 PM
System		File Fold...	1/1/06 2:00 PM
TitanCals		File Fold...	10/27/14 12:26 PM
BrukerInstaller.xls	1 KB	Microsof...	10/21/14 11:35 AM
BrukerInstallerLog0.txt	141 KB	Text Do...	12/6/16 3:04 AM
Cal DualMining Context ...	8 KB	SNPSH...	5/...
Cargill.aen	5 KB	AEN File	6/...
eventlog0.txt	1,505 KB	Text Do...	1/...
eventlog1.txt	721 KB	Text Do...	1/...
eventlog2.txt	931 KB	Text Do...	1/...
GeoChem.aen	22 KB	AEN File	4/...
InstrumentDef.idf	34 KB	IDF File	1/...
MudrockMajor.aen	5 KB	AEN File	7/...
Mudrock-Majors.aen	5 KB	AEN File	5/...
MudrockTrace.aen	5 KB	AEN File	7/...
Mudrock-Trace.aen	5 KB	AEN File	5/...
Obsidian.aen	5 KB	AEN File	7/...
plantCal.aen	5 KB	AEN File	12/...
PreciousMetals.aen	19 KB	AEN File	11/...
RoHS.alt	47 KB	ALT File	11/...
SmartGrade.aen	52 KB	AEN File	4/...
Soil.aen	21 KB	AEN File	4/...
SOILTEST50.aen	5 KB	AEN File	4/...
StdAlloy.aen	1,411 KB	AEN File	4/...
tracelog0.txt	523 KB	Text Do...	1/1/06 9:01 PM

Browse For Folder

- Desktop
  - Libraries
  - Homegroup
  - lee
  - Computer
  - Network
  - Control Panel
  - Recycle Bin
  - 3-26-17\_STDS
  - 4.2 ky event
  - Si DUGWAY OBS calibration

Make New Folder OK Cancel

In the window that opens, navigate to where you would like to download the data

- My Computer
  - PRL (A:)
  - (C:)
  - \\Mac\Google Drive\ (V:)
  - \\Mac\Dropbox\ (W:)
  - \\Mac\Photo Library\ (X:)
  - \\Mac\iCloud\ (Y:)
  - \\Mac\Home\ (Z:)



\BRUKER

- 800N2682
  - Application Data
  - ARCHIVE
  - BINARIES
  - BRUKER
    - Autold
    - Data
    - Documents and Settings
    - EDX Databases
    - lost+found
    - package
    - Setup
    - System
    - TitanCals
  - LBOOT
  - My Documents
  - Network
  - Program Files
  - Recycled
  - Temp

Name	Size	Type	Date Modified
Autold		File Fold...	1/1/06 2:00 PM
Data		File Fold...	10/21/14 11:35 AM
Documents and Settings		File Fold...	1/1/06 6:00 AM
EDX Databases		File Fold...	1/1/06 2:00 PM
lost+found		File Fold...	1/11/06 12:59 PM
package		File Fold...	1/1/06 2:00 PM
Setup		File Fold...	1/1/06 2:00 PM
System		File Fold...	1/1/06 2:00 PM
TitanCals		File Fold...	10/27/14 12:26 PM
BrukerInstaller.xls	1 KB	Microsof...	10/21/14 11:35 AM
BrukerInstallerLog0.txt	141 KB	Text Do...	12/6/16 3:04 AM
Cal DualMining Context ...	8 KB	SNPSH...	5/30/16 8:10 AM
Cargill.aen	5 KB	AEN File	6/20/16 11:08 AM
eventlog0.txt	1,505 KB	Text Do...	
eventlog1.txt	721 KB	Text Do...	
eventlog2.txt	931 KB	Text Do...	
GeoChem.aen	22 KB	AEN File	
InstrumentDef.idf	34 KB	IDF File	
MudrockMajor.aen	5 KB	AEN File	
Mudrock-Majors.aen	5 KB	AEN File	
MudrockTrace.aen	5 KB	AEN File	
Mudrock-Trace.aen	5 KB	AEN File	
Obsidian.aen	5 KB	AEN File	
plantCal.aen	5 KB	AEN File	
PreciousMetals.aen	19 KB	AEN File	
RoHS.alt	47 KB	ALT File	11/2/15 2:04 PM
SmartGrade.aen	52 KB	AEN File	4/11/16 10:21 AM
Soil.aen	21 KB	AEN File	4/11/16 10:21 AM
SOILTEST50.aen	5 KB	AEN File	4/30/16 4:18 PM
StdAlloy.aen	1,411 KB	AEN File	4/11/16 10:22 AM
tracelog0.txt	523 KB	Text Do...	1/1/06 9:01 PM

Wait until the data is done copying, and you are good to go!

Downloading...

\BRUKER\Data\00083-Spectrum Only.pdz -> \\Mac\Home\...

( 61/1456 ) : 8788 of 8788 bytes copied.

Cancel

- My Computer
  - PRL (A:)
  - (C:)
  - \\Mac\Google Drive\ (V:)
  - \\Mac\Dropbox\ (W:)
  - \\Mac\Photo Library\ (X:)
  - \\Mac\iCloud\ (Y:)
  - \\Mac\Home\ (Z:)