FILTER VOLTAGE AND CURRENT FOR XRF SPECTROSCOPY USER GUIDE FOR ANALYZING COMMON ELEMENTS PRESENT IN PHOTOGRAPH AND PAPER COLLECTIONS

XRF IDENTIFICATION OF SILVER

Element: Ag

- 1. Insert green filter
- 2. 45 kV
- 3. 12 micro amps
- 4. No vacuum

XRF IDENTIFICATION OF PLATINUM

Element: Pt

- 1. Insert green filter
- 2. 45 kV
- 3. 12 micro amps
- 4. No vacuum

XRF IDENTIFICATION OF GOLD

Element: Ag

Useful when analyzing photographs to observe presence of gold toning

- 1. Insert green filter
- 2. 40 kV
- 3. 12 micro amps
- 4. No vacuum

XRF IDENTIFICATION OF PALLADIUM

Element: Pd

- 1. Insert green filter
- 2. 45 kV
- 3. 12 micro amps
- 4. No vacuum

XRF IDENTIFICATION OF IRON

Element: Fe

Useful when analyzing cyanotypes and toned gelatin silver photographs (exhibit bluish tint)

- 1. Insert yellow filter
- 2. 40 kV
- 3. 1.0 4.0 micro amps
- 4. No vacuum

XRF IDENTIFICATION OF SULFUR

Element: S

May be useful when comparing albumen photographs to lightly albumenized salted paper prints. May be used to indicate sulfur-toned gelatin silver prints.

- 1. Insert blue filter
- 2. 15 or 20 kV
- 3. 15 or higher micro amps
- 4. Use vacuum
- 5. Attach helium cap

XRF IDENTIFICATION OF BARIUM

Element: Ba

Useful when analyzing photographs to observe presence of baryta layer; may be useful when calculating barium/strontium ratio to determine source of barium to help date photographs

- 1. No filter
- 2. 15 kV
- 3. 15 micro amps
- 4. Vacuum

XRF IDENTIFICATION OF LEAD

Element: Pb

Common additive to adhesives; lead white paint

- 1. Insert red filter
- 2. 40 kV
- 3. 4.0 to 8.0 micro amps
- 4. No vacuum

XRF IDENTIFICATION OF MERCURY

Element: Hg

- 1. Insert green filter
- 2. 40 kV
- 3. 12 micro amps
- 4. No vacuum

XRF IDENTIFICATION OF SELENIUM

Element: Se

Useful when analyzing photographs to observe presence of selenium toning

- 1. Insert yellow filter
- 2. 40 kV
- 3. 1.0 to 4.0 micro amps
- 4. No vacuum

XRF IDENTIFICATION OF TITANIUM

Element: Ti

- 1. Insert yellow filter
- 2. 40 kV
- 3. 1.0 to 4.0 micro amps
- 4. No vacuum

XRF IDENTIFICATION OF URANIUM

Useful when analyzing yellow toned photographs for presence of uranium toning $Element: \ U$

- 1. Insert red filter
- 2. 40 kV
- 3. 4.0 to 8.0 micro amps
- 4. No vacuum

XRF IDENTIFICATION OF COPPER

Useful when analyzing reddish-tinted photographs for presence of copper toning

Element: Cu

- 1. Insert yellow filter
- 2. 40 kV
- 3. 1.0 to 4.0 micro amps
- 4. No vacuum