



Ready to Test

APPLICATION  
plantCal

SETTINGS

METHOD  
plantCal

DISPLAY  
(None)

Camera

Utilities

Edit Info

Logoff

Results

Backup Data

Library  
Maintenance

System Setup

Remote Control

File Explorer

Back



28



Applications

Camera Setup

Connections

Date\Time

Languages

Output Control

Run Orders

Safety

More

Back



29



### Connection

- Enable Wireless
- Enable Bluetooth
- Enable USB
- Enable Serial

Configure Connections

OK

Back

[Bluetooth] [USB] [Wi-Fi]

Enable

Instrument Name

Scan for Devices

Device	Paired	Se'l'd

OK

Back



Enable  IP: 192.168.0.21

Network Name (SSID)	P	C
Add New...		
NETGEAR21	N	N
La Casa de Biologia	Y	Y
Casa De Paco	N	N
xfinitywifi	N	N

Connected to La Casa de Biologia

Connect IP Setup Delete

OK

Back

File Trigger Unknown Display Help

31

## Wireless Network Properties

Network Name (SSID)

Set as Adhoc Network Connection

Wireless network key (WEP)

This network requires a key for:

Authentication:

Encryption:

Network key:

Key index:

The key is provided automatically

IEEE 802.1X Authentication

Enable 802.1X authentication

EAP Type:

File Trigger Unknown Display Help

31

## Wireless Network Properties

Network Name (SSID)

Set as Adhoc Network Connection

Wireless network key (WEP)

This network requires a key for:

Authentication:

Encryption:

Network key:

Key index:

The key is protected

IEEE 802.1X Authentication

Enable 802.1X authentication

EAP Type:

OK Back





(8)

31



# Wireless Network Properties

Network Name (SSID)

Set as Adhoc Network Connection

## Wireless network key (WEP)

This network requires a key for:

Authentication:

Encryption:

Network key:

Key index:

The key is protected

The key is protected

The key is protected

## IEEE 802.1X Authentication

Enable 802.1X authentication

EAP Type:

OK

Back

File Trigger Unknown Display Help

32

## Wireless Network Properties

Network Name (SSID)

Set as Adhoc Network Connection

Wireless network key (WEP)

This network requires a key for:

Authentication:

Encryption:

Network key:

Key index:

The key is provided automatically

IEEE 802.1X Authentication

Enable 802.1X authentication

EAP Type:

OK Back

File Trigger Unknown Display Help

32

## Wireless Network Properties

Network Name (SSID)

Set as Adhoc Network Connection

Wireless network key (WEP)

This network requires a key for:

Authentication:

Encryption:

Network key:

Key index:

The key is provided automatically

IEEE 802.1X Authentication

Enable 802.1X authentication

EAP Type:

OK Back



# Network Key

YourPassword|

1	2	3	4	5	6	7	8	9	0
q	w	e	r	t	y	u	i	o	p
a	s	d	f	g	h	j	k	l	;
CAPS	z	x	c	v	b	n	m	Shift	
	-	=	[	]	\		.	/	
Delete		Clear		Backspace			Reset		
OK		LT	Space		RT	Cancel			

File Trigger Unknown Display Help

32

## Wireless Network Properties

Network Name (SSID)

Set as Adhoc Network Connection

Wireless network key (WEP)

This network requires a key for:

Authentication:

Encryption:

Network key:

Key index:

The key is provided automatically

IEEE 802.1X Authentication

Enable 802.1X authentication

EAP Type:

OK Back



Enable  IP: 192.168.0.21

Network Name (SSID)	P	C
Add New...		
NETGEAR21	N	N
La Casa de Biologia	Y	Y
Casa De Paco	N	N
xfinitywifi	N	N

Connected to La Casa de Biologia

Connect IP Setup Delete

OK Back



33



### WiFi IP Setup

DHCP

IP 192 168 0 21

Mask 255 255 255 0

Gateway 192 168 0 1

MAC 00:08:EE:0A:C2:E2

Renew

OK

Back



33



### Connection

- Enable Wireless
- Enable Bluetooth
- Enable USB
- Enable Serial

Configure Connections

OK

Back



Password (None) X

User: test ▼

Password:

Change Password Ok Cancel